

96-Well Magnet Plate LE400

Low Elution PCR



Catalog # (P/N): LE400

Precautions:

CAUTION - STRONG MAGNETIC FIELD. This Permagen Labware 96-well Magnet plate contains very strong neodymium iron boron (NdFeB) magnets. People with pacemakers or other forms of Implantable cardioverter-defibrillators (ICD's) should always avoid this product as it may interfere with its operation. Never allow this product to come into contact with other magnets, metal, magnetized media such as credit cards etc., as damage and or injury can occur. Do not Autoclave or heat above 80°C, doing so will damage the magnets.

USER INFORMATION

Product Description:

Permagen Labware's 96-well Magnet Plate is designed for manual or automated paramagnetic bead separation with many standard 96-well PCR plates. Flat PCR plates are recommended for optimal performance. The magnets in our LE400 Magnet plate will form beads to the side of each well leaving a clean area for aspiration. Separate as low as 5 µL. Standard SBS footprint base which will fit into almost any standard liquid handler.

Care/ Cleaning:

If solutions are spilled onto the magnet plate carefully rinse with tap water and allow to air dry completely before use. A mild soap may also be used if more extreme cleaning is necessary. Always use care when storing product when not in use.

OTHER INFORMATION

Warranty and Liability:

For research use only. Not for use in diagnostic procedures.

Permagen Labware, LTD. Is focused on delivering industry leading product quality, and performance. Please visit www.permagenlabware.com for full warranty terms and conditions.

Any information contained herein is subject to change without notice. Permagen Labware, LTD. assumes no responsibility for any errors that may appear in this document.

Permagen Labware is a registered trademark of Permagen Labware, LTD
© 2015 Permagen Labware, LTD. All rights reserved.